

Product Information

Peroxisome Isolation Kit

Product Code **PEROX1**

Storage Temperature 2–8 °C

TECHNICAL BULLETIN

Product Description

Peroxisomes are organelles ubiquitously distributed in most prokaryotic and eukaryotic cells. They are spherical particles with a diameter of 0.2–1.0 μm . Peroxisomes are subcompartmentalized with a unit membrane surrounding a matrix compartment, similar to mitochondria, that contains the bulk of the peroxisomal enzymes. Peroxisomes are most abundant in the mammalian liver and kidney.^{1–3} In hepatocytes and kidney cortex epithelial cells, peroxisomes are large (0.3–1.5 μm); whereas, in other cells such as fibroblasts or smooth muscle cells, the peroxisomes are small (0.05–0.2 μm). Isolated peroxisomes are used for studying lipid β -oxidation,⁴ amino acid metabolism,⁵ and biosynthesis of ether-linked glycerolipids⁶ and bile acids.^{7,8}

The Peroxisome Isolation Kit provides all necessary reagents and a comprehensive protocol for isolating peroxisomes from animal tissues and from cultured cells by differential centrifugation followed by density gradient centrifugation. The kit was tested with rat liver and kidney, rabbit liver, and with HEK293 and HepG2 cell lines.

The presence of peroxisomes can be determined by measuring the activity of the peroxisomal enzyme marker, catalase (Catalase Assay Kit, Product Code CAT100). In addition, the detection of peroxisomal membrane proteins and their degree of enrichment in the purified fraction can be performed by immunoblot, using antibodies specific for the peroxisomal membrane specific marker PMP70 (Product Code P 0497). Separation of peroxisomes from mitochondria can be determined by measuring the activity of the mitochondrial enzyme marker, cytochrome c oxidase (Cytochrome c Oxidase Assay Kit, Product Code CYTOCOX1).

Components

The kit is sufficient for the preparation of peroxisomes from 50 g of tissue or ~20 ml of packed cells

Peroxisome Extraction Buffer 5x 100 ml
Product Code P 7247
25 mM MOPS, pH 7.65, with 1.25 M sucrose,
5 mM EDTA, and 0.5% ethanol

Optiprep Dilution Buffer 20x 10 ml
Product Code O 4889
100 mM MOPS, pH 8.0, with 20 mM EDTA and
2% ethanol

Protease Inhibitor Cocktail for use with 5 ml
mammalian cell and tissue extracts
Product Code P 8340

OptiPrep Density Gradient Medium 100 ml
Product Code D 1556
60% (w/v) solution of iodixanol in water

Equipment and Reagents Required But Not Provided

- Sorvall[®] RC-5C centrifuge with SS-34 head or equivalent
- Ultracentrifuge and a fixed angle head suitable for centrifugation at 100,000 x g with 8 ml sample volume per tube
- Ultrapure water
- Dulbecco's Phosphate Buffered Saline (PBS) (Product Code D 8537)
- Microcentrifuge tubes
- Pasteur pipettes

Equipment Required for Peroxisome Preparation from Tissue Extracts

- Homogenizers:
 - For initial homogenization – Tissue homogenizer (Ultra-Turrax® T-25 with S25N 18G head or equivalent)
 - For homogenization of 25,000 x g pellet - Overhead electric motor together with Potter-Elvehjem PTFE pestle in glass tube homogenizer - 3 ml (Product Code P 7734) or 8 ml (Product Code P 7859)
- Scalpel, forceps, and glass plate
- 40 ml polypropylene centrifuge tubes

Equipment and Reagents Required for Peroxisome Preparation from Cell Culture Extracts

- Trypsin-EDTA (for removal of adherent cells from vessel surface, Product Code T 4049)
- Trypan Blue (Product Code T 8154)
- Sorvall RT-6000B centrifuge or equivalent
- Adaptor for microcentrifuge tubes in Sorvall RC-5C
- 14 ml centrifuge tubes with adaptor for SS-34 head
- Homogenizers:
 - For initial homogenization - 7 ml Dounce glass tissue grinder (Product Code T 0566) with small clearance pestle (Product Code P 1235)
 - For homogenization of 25,000 x g pellet: Pellet pestle (Product Code Z35,994-7) and motor for pellet pestle (Product Code Z35,997-1).

Additional Equipment and Reagents Required to Measure Enzyme Activities

- Spectrophotometer with a 1 ml cuvette (for measurement of enzyme activities)
- Catalase Assay Kit (Product Code CAT100)
- Cytochrome c Oxidase Assay Kit (Product Code CYTOCOX1).

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

It is recommended to use ultrapure (17 MΩ-cm or equivalent) water when preparing the reagents.

1x Peroxisome Extraction Buffer - Dilute an aliquot of the Peroxisome Extraction Buffer 5x (Product Code P 7247) 5-fold with ultrapure water. Keep the diluted Peroxisome Extraction Buffer at 4 °C until use. Just before use add Protease Inhibitor Cocktail for mammalian cell and tissue extracts (Product Code P 8340) to the diluted Peroxisome Extraction Buffer at a concentration of 1% (v/v). The diluted Peroxisome Extraction Buffer with the 1% (v/v) Protease Inhibitor Cocktail is the 1x Peroxisome Extraction Buffer.

Suggested volumes of 1x Peroxisome Extraction Buffer:

- For tissue extracts: use a minimal tissue weight of 4 g and prepare 25 ml of buffer.
- For cell culture extracts: prepare a minimum of 2×10^8 cells and 10 ml of buffer.

1x Optiprep Dilution Buffer - Dilute an aliquot of the Optiprep Dilution Buffer 20x (Product Code O 4889) 20-fold with water. Keep the 1x Optiprep Dilution Buffer at 4 °C until use. Suggested volumes of 1x Optiprep Dilution Buffer:

- For tissue extracts from 4 g tissue: 25 ml.
- For cell culture extracts of at least 2×10^8 cells: 20 ml.

OptiPrep Density Gradient Medium Solutions - For separation in 8 ml centrifuge tubes prepare:

- 27.5% Optiprep Solution – Mix 4.58 ml of OptiPrep Density Gradient Medium [60% (w/v)] with 5.42 ml of 1x Optiprep Dilution Buffer. Mix well by inversion.
- 20% Optiprep Solution – Mix 3.33 ml of OptiPrep Density Gradient Medium [60% (w/v)] with 6.67 ml of 1x Optiprep Dilution Buffer. Mix well by inversion.

Storage/Stability

The kit is shipped on wet ice and storage at 2–8 °C is recommended. Upon receiving the kit, the Protease Inhibitor Cocktail (Product Code P 8340) should be stored at –20 °C and the OptiPrep Density Gradient Medium (Product Code D 1556) should be stored at room temperature. When stored properly, the components in this kit are stable for 24 months.

Procedure

Peroxisomes with different degrees of purity can be easily prepared from animal tissues using a simple method of homogenization with the aid of an Ultra-Turrax T25 tissue homogenizer followed by differential centrifugation. Tissue culture cells need to be homogenized in a Dounce homogenizer and the degree of breakage carefully followed by Trypan Blue staining. Excess breakage will cause widespread damage to the peroxisomes.

The serial centrifugations include:

- Low speed centrifugation (1000 x g)
- Low speed centrifugation (2000 x g)
- Medium speed centrifugation (25,000 x g)

The serial centrifugations remove nuclei, cell debris, mitochondria, and lipids to obtain a crude peroxisomal fraction (CPF). The CPF pellet is the starting material for the preparation of purified peroxisomes.

A highly enriched peroxisome fraction can be prepared from the CPF pellet by efficient removal of the other organelles using an Optiprep density gradient: the CPF suspended in 22.5% Optiprep is layered between 27.5% and 20% Optiprep layers, and centrifuged at 100,00 x g for 1.5 hours. The purified peroxisome fraction is collected from the bottom of the tube.⁹

Note: If an ultracentrifuge is not available it is possible to centrifuge the same tubes in a Sorvall RC-5C centrifuge for 3.5 hours at 45,000 x g (19,400 rpm in an SS-34 head).

A flow diagram for the various peroxisome preparations is shown in the Appendix.

A. Preparation of Crude Peroxisomal Fraction (CPF) from Animal Tissue (~4 grams of tissue)

Perform the procedure at 4 °C. All solutions and equipment should be cooled before use. Homogenize the samples using an UltraTurrax T-25 tissue homogenizer with S25N 18G head.

1. Use a fresh tissue sample from an animal that was starved overnight and sacrificed the next morning.
2. Wash the tissue sample three times with 10-15 ml of ice cold PBS by placing the tissue in a dish, shaking gently for few minutes, and removing the PBS. Place the tissue on a paper towel in order to absorb excess liquid and blood clots, if present. Cut the tissue into small pieces (1.5 - 2 cm) and repeat the wash step.

Note: This step is performed in order to remove excess blood from certain tissues (e.g. liver).

3. Blot the tissue on a paper towel and weigh.
4. Cut the tissue on a glass plate with the aid of a scalpel into small slices (0.3-0.5 cm in width). Transfer the slices into a 40 ml polypropylene centrifuge tube. Add 4 volumes of the 1x Peroxisome Extraction Buffer per gram of tissue (i.e. 16 ml per 4 grams), and homogenize the sample as follows: Homogenization at 8,000 rpm for 5 seconds followed by 2 additional homogenizations at 9,500 rpm for 5 seconds each.
5. Wash the homogenizer head with 1 ml of the 1x Peroxisome Extraction Buffer and add the wash to the previous homogenate. Keep the homogenate on ice.
6. Centrifuge the homogenate at 1,000 x g for 10 minutes at 4 °C. Remove the floating lipid layer by aspiration, transfer the supernatant to another centrifuge tube using a pipette, and discard the pellet. This step removes nuclei and other cell debris.
Note: To monitor the degree of purification following the different centrifugation steps, it is recommended to save a sample (100-200 µl) of the 1,000 x g supernatant for subsequent assays.
7. Centrifuge at 2,000 x g for 10 minutes at 4 °C. Remove the floating lipid layer by aspiration, transfer the supernatant to another centrifuge tube using a pipette, and discard the pellet. This step removes heavy mitochondria.
8. Centrifuge at 25,000 x g for 20 minutes at 4 °C. Aspirate off the supernatant liquid and resuspend the pellet in a minimal volume of 1x Peroxisome Extraction Buffer. It is recommended to use 0.4 ml per gram of original tissue (i.e. 1.6 ml per 4 grams).

The suspension obtained in step 8 is the Crude Peroxisomal Fraction (**CPF**) and contains a mixture of light mitochondria, lysosomes, peroxisomes, and endoplasmic reticulum (ER).

Note: The CPF may be kept overnight at 4 °C and then separated the next day on a density gradient.

B. Preparation of Crude Peroxisomal Fraction (CPF) from cell cultures (~2 x 10⁸ cells)

Perform the procedure at 4 °C. All solutions and equipment should be cooled before use. Homogenize the samples using a 7 ml Dounce glass tissue grinder (Product Code T 0566) with small clearance pestle (Product Code P 1235).

Note: This procedure requires a relatively large amount of cells, 1-2 ml packed cell volume representing at least 2 x 10⁸ cells.

1. Wash the cells

For adherent cells:

- Wash cells with PBS to remove the serum (~8 ml per 100 cm²).
- Detach the cells from the surface with Trypsin or Trypsin-EDTA.
- Pour the trypsinized cells into growth medium containing 10% Fetal Bovine Serum (Product Code F 6178) (4 ml per 100 cm²). This quenches the proteolytic action of the trypsin.

For cells in suspension, start with step 2.

2. Centrifuge the cell suspension for 5 minutes at 250 x g (~1,200 rpm in a Sorvall RT-6000B centrifuge). Discard the supernatant and resuspend the cells in 40 ml of PBS.
3. Centrifuge the resuspended cells for 5 minutes at 250 x g and discard the supernatant. Wash the cells twice in this fashion. The packed cell volume (PCV) should be 1.5-3 ml.
4. Add 2.7 PCV of 1x Peroxisome Extraction Buffer and vortex to achieve an even suspension.
5. Break the cells in a 7 ml Dounce homogenizer (Product Code T 0566) using Pestle B (small clearance). This may necessitate splitting the fraction into two portions.
6. After every 5 strokes with the pestle, stain an aliquot of cells with Trypan Blue and examine under a microscope to assess the degree of breakage. Normally 15-25 strokes are sufficient to achieve 80-85% breakage. Do not try to achieve higher breakage levels as this will lead to severe damage of the peroxisomes.
7. Centrifuge the sample at 1,000 x g for 10 minutes.
8. Transfer the supernatant to a new centrifuge tube.
Note: To monitor the degree of purification following the different centrifugation steps, it is recommended to save a sample (100-200 µl) of the 1,000 x g supernatant for subsequent assays.
9. Centrifuge the sample at 2,000 x g for 10 minutes.
10. Transfer the supernatant to a new centrifuge tube.
11. Centrifuge the sample at 25,000 x g for 20 minutes in Eppendorf® tubes using a special adaptor for the Sorvall centrifuge.
12. Remove the supernatant liquid and collect the pellet in a minimal volume of 1x Peroxisome Extraction Buffer (~0.4 ml per 10⁸ cells).
13. Uniformly suspend the pellet in a single Eppendorf tube by using a pellet pestle.

The suspension obtained in step 13 is the Crude Peroxisomal Fraction (**CPF**) and contains a mixture of light mitochondria, lysosomes, peroxisomes, and endoplasmic reticulum.

Note: The CPF may be kept overnight at 4 °C and then separated the next day on a density gradient.

C. Isolation of Peroxisomes on a Density Gradient

The objective of density gradient centrifugation is to obtain enriched peroxisomes separated from other organelles, e.g. mitochondria.

This procedure uses an 8 ml ultracentrifuge tube.

1. Dilute the CPF (section A, step 8 or section B, step 13) as follows: to 1.2 ml of the CPF (crude peroxisomal fraction) add 1.69 ml of the OptiPrep Density Gradient Medium and 1.61 ml of the 1x Optiprep Dilution Buffer. This yields a diluted CPF sample with a final volume of 4.5 ml and an OptiPrep concentration of 22.5%. If the CPF sample is not properly diluted, it will tend to drop into the lower density layer.
2. Place 2 ml of the 27.5% Optiprep Solution into an 8 ml ultracentrifuge tube using a Pasteur pipette and then overlay with 4.0 ml of the diluted CPF sample. Overlay the sample with 2 ml of the 20% Optiprep Solution.
Note: Overlay the different layers by carefully applying the solutions onto the wall of the tube.
3. Close the tube and ensure that it is balanced against another tube.
4. Centrifuge for 1.5 hours at 100,000 x g. When deaccelerating the centrifuge, brake to 800 rpm and then let it to come to a stop without use of the brake. The tube will show a cloudy, floating layer on the top of the gradient that mainly consists of ER and lysosomes, a ring at the 22.5%/27.5% interface that is mitochondria, and a slightly floating pellet in the bottom layer.
5. Carefully aspirate off the top layers including the ring at the 22.5%/27.5% interface. This step is important in order to obtain a fraction that is as free as possible of mitochondria.
6. Withdraw the bottom layer containing the **purified peroxisomes**.

Note: The purified peroxisomes may also be kept at 4 °C for up to 24 hours with little degradation. If an immunoblot of specific proteins in the peroxisome is desired, it is advisable to add the sample buffer immediately after preparation of the purified fraction and then to freeze the sample.

D. Analysis of Peroxisome Purification

The degree of purification of the peroxisomes can be evaluated by comparing the protein concentration, catalase (peroxisomal marker) and cytochrome c oxidase (mitochondrial marker) activities of the purified peroxisomes to that of the 1000 x *g* supernatant. The ratio of catalase activity to cytochrome c oxidase activity in the purified peroxisomes should be at least 15-30 times greater than that found in the 1,000 x *g* supernatant.

Catalase activity must be determined by a colorimetric method since Optiprep interferes with the UV method.

The purified peroxisomes can be assayed according to the following guidelines:

- a. Protein concentration by Bradford method:
Dilute 10-fold and use 10-40 µl per test.
- b. Catalase activity:
For tissues, dilute 20 to 100-fold and use 2.5-5.0 µl per test.
For cell extracts, dilute 4-fold and use 2.5-5.0 µl per test.
- c. Cytochrome c oxidase activity:
Dilute 3-fold and use 20-100 µl per test.

The 1000 x *g* supernatant can be assayed according to the following guidelines:


- a. For protein determination: Dilute 100-fold and use 20-40 µl per test.
- b. Catalase activity:
For tissues, dilute 100-fold and use 2.5-5.0 µl per test.
For cell extracts, dilute 6-fold and use 2.5-5.0 µl per test.
- c. Cytochrome c oxidase activity: Dilute 3-fold and use 5-10 µl per test.

Results

An immunoblot of the purified fractions using the anti-PMP70 antibody (Product Code P0497), an antibody against peroxisomal membrane protein 70 (PMP70), is shown below in Figure 1.

Figure 1.

Enrichment of peroxisomes in isolated fractions detected by anti-PMP70 antibody

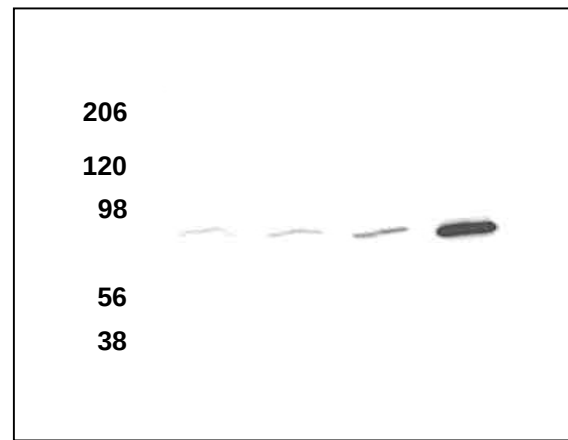

Rat liver peroxisomes were purified according to the isolation procedure. Samples of 0.2 µg protein from the different purification stages/fractions were run on 7.5% SDS-PAGE. Immunoblotting was performed using anti-PMP70 at a dilution of 1:2000.

Lane A: 1,000 x *g* supernatant

Lane B: The endoplasmic reticulum fraction from the Optiprep gradient

Lane C: The fraction containing mainly mitochondria (with some peroxisomes) from the Optiprep gradient

Lane D: The purified peroxisome fraction from the Optiprep gradient.


References

1. Hartl, F-U., *et al.*, Improved isolation and purification of rat liver peroxisomes by combined rate zonal and equilibrium density centrifugation. *Arch. Biochem. Biophys.*, **237**, 124-134 (1985).
2. Volkl, A., and Fahimi, H.D., Isolation and characterization of peroxisomes from the liver of normal untreated rats. *Eur. J. Biochem.*, **149**, 257-265 (1985).
3. Zaar, K., Structure and function of peroxisomes in the mammalian kidney. *Eur. J. Cell Biol.*, **59**, 233-254 (1992).
4. Lazarow, P.B., and De Duve, C., A fatty acyl-CoA oxidizing system in rat liver peroxisomes; enhancement by clofibrate a hypolipidemic drug. *Proc. Natl. Acad. Sci. USA*, **73**, 2043-2046 (1976).
5. Mannaerts, G.P., and Van Veldhoven, P.P., Metabolic pathways in mammalian peroxisomes. *Biochimie*, **75**, 147-158 (1993).
6. Burdett, K., *et al.*, Peroxisomal localization of Acyl-coenzyme A reductase (long chain alcohol forming) in guinea pig intestine mucosal cells. *J. Biol. Chem.*, **266**, 12201-12206 (1991).
7. Rodrigues, C.M.P., *et al.*, Formation of Δ^{22} -bile acids in rats is not gender specific and occurs in the peroxisomes. *J. Lipid Res.*, **37**, 540-550 (1996).
8. Une, M., *et al.*, Comparison of side chain oxidation of potential C₂₇-bile acid intermediates between mitochondria and peroxisomes of the rat liver: presence of β -oxidation activity for bile acid biosynthesis in mitochondria. *J. Lipid Res.*, **37**, 2550-2556 (1996).
9. Graham, J., *et al.*, The preparation of subcellular organelles from mouse liver in self-generated gradients of Iodixanol. *Anal. Biochem.*, **220**, 367-373 (1994).

Sorvall is a registered trademark of Kendro Laboratory Products.

Ultra-Turrax is a registered trademark of IKA WORKS, Inc.

Eppendorf is a registered trademark of Eppendorf-Netheler-Hinz GmbH.

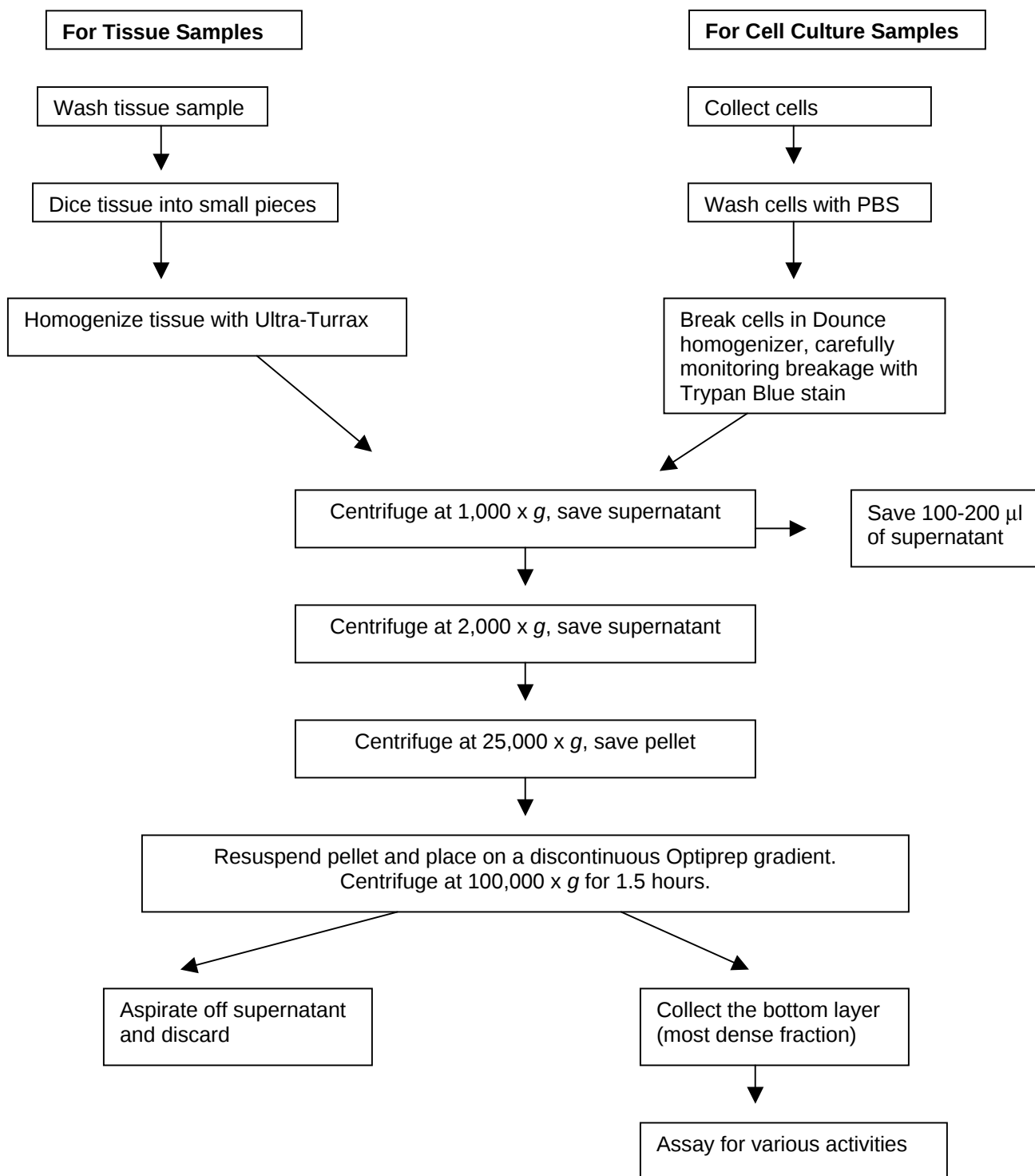Appendix – Flow Chart of Peroxisome Preparation

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.